

Mixing Chamber Assembly Instructions:


Parts List:

- 1) Chamber
- 2) Input port (smaller dia, near back of chamber)
- 3) Output port (larger dia, near the lid)
- 4) Metal washers (4)
- 5) Plastic Washers (4)
- 6) Nut (2)
- 7) Luer Connectors (2).

1) Screw the Luer Connectors to the top of the chamber.



1: Luer Connectors

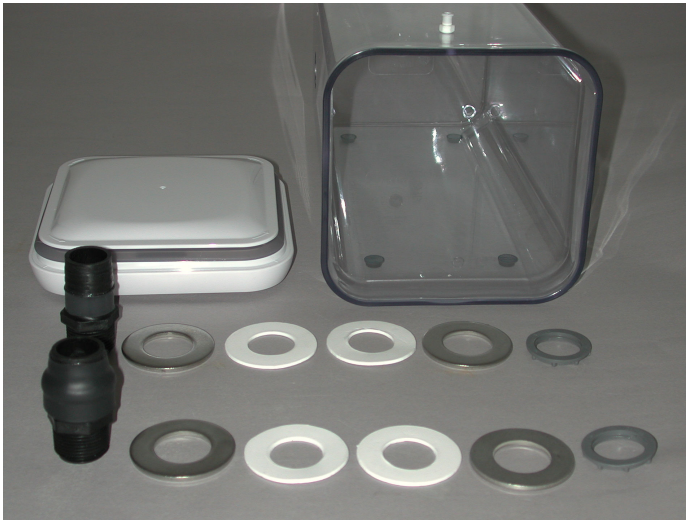


2: parts



3: Metal and Plastic washer on port

2) Place the metal washer and then a plastic washer on each port. Insert it through the hole in the chamber wall and place the plastic washer, then the metal washer and then the nut, from the inside. Hand Tighten, both ports.

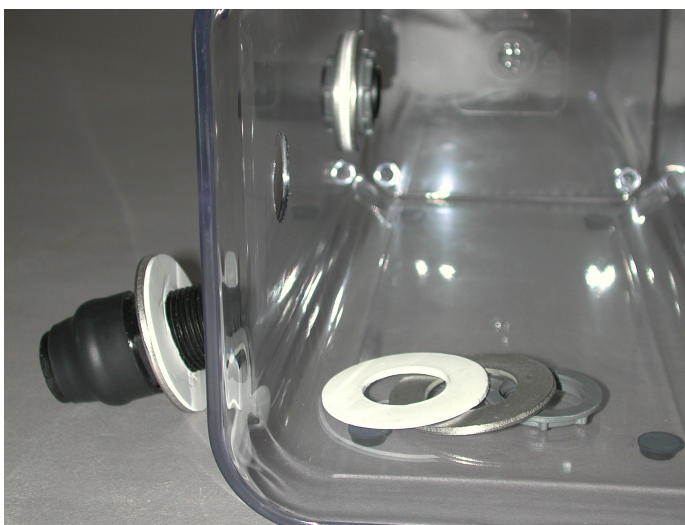

4: input port assembled and output port partially assembled



5: Assembled Chamber.

3) The chamber is now ready for use.